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Nr./Art. 85754

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Instructions Stethoscope

Nr./Art. 85754

Important – Please read instructions carefully and keep for future reference!

Using the stethoscope correctly:

- ▶ The stethoscope is intended exclusively for educational use.
- ▶ The stethoscope is not intended for medical purposes.
- ▶ The stethoscope is not a toy.
- ▶ Children should only use the stethoscope under adult supervision.
- ▶ The stethoscope has small parts that could be swallowed, and is therefore not suitable for children under the age of three years.
- ▶ Always clean the earpieces with a damp cloth before and after use.

Warnings:

- ▶ Before use, check that all parts are in perfect condition. Never use damaged parts.
- ▶ The stethoscope is not a toy. It is not suitable for children under the age of three years.

Parts and material:

The stethoscope consists of three parts.

The binaurals with the earpieces consist of seamless brass tubes.

The high-grade steel spring which is fastened to the frame ensures suitable elasticity and a perfect fit.

The smooth plastic earpieces sit comfortably in the ear.

The thick tubing is made of PVC.

The chestpiece is also made of brass.

All the stethoscope's components are free of nickel and latex.

Storage:

- ▶ Keep the stethoscope out of reach of children.
- ▶ Keep the stethoscope in the supplied box.
- ▶ Keep the stethoscope in a dry place and protect it against the effects of direct sunlight.
- ▶ Do not keep the stethoscope near heat sources.

Preparations for use:

Push the two ends of the Y-shaped tubing as far as possible onto the metal ends of the binaurals.

Push the other end of the tubing as far as possible onto the stem of the chestpiece.

Using the stethoscope correctly and safely:

The two earpieces should be placed loosely in the ear, blocking the auditory canal to prevent distraction from external sounds.

Make sure that the earpiece does not enter the auditory canal, as it could damage it.

The chestpiece can now be held against the area of the body to which you wish to listen.

How the stethoscope works:

The stethoscope's chestpiece contains a membrane that picks up acoustic vibrations and starts vibrating itself.

The membrane conducts these vibrations to the air column in the stethoscope's tubing.

The vibrations are conveyed via the binaurals to the user's eardrums.

The user thus hears the sounds in amplified form.

Maintenance and cleaning:

After use, clean the components with a damp cloth only. Do not use detergent. Make sure that all components are completely dry before they are used again. Always clean the earpieces with a damp cloth before and after use.

What to do with the packaging:

Make sure that the packaging is disposed of immediately after removal in a way that does not harm the environment.

The plastic bag presents a suffocation hazard for young children.

Warranty:

The legal warranty period of two years applies to this product.

This period starts on the date of product delivery.

Please contact directly your supplier for any warranty-related matters.

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Arnulf Betzold GmbH
Betzold Lernmedien GmbH